Datasheet

HSA-X-S-1G4-SI

Ultra High Speed Photoreceiver with Si-PIN Photodiode



The picture shows model HSA-X-S-1G4-SI-FST


Features

- Si-PIN photodiode
- Bandwidth 10 kHz 1.4 GHz
- Amplifier transimpedance gain 5.0 × 10³ V/A
- Max. conversion gain 2.55 × 10³ V/W @ 760 nm
- Spectral range 320 1000 nm
- Free-space input 1.035"-40 threaded, easily convertible to fiber optic input (FC and FSMA) with optionally available screw-on adapters
- Fiber optic input also available as permanently mounted FC-input
- UNC 8-32 and M4 tapped holes for mounting on standard posts with metric and imperial thread

Applications

- Spectroscopy
- Ultra-fast pulse and transient measurements
- Optical triggering
- Optical front-end for oscilloscopes and ultra-fast A/D converters

Block Diagram

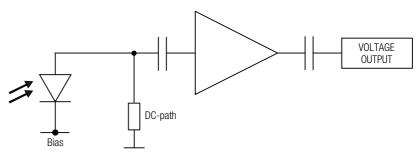

BS01-HSA-X-S_R01

Intended Use

The HSA-X-S-1G4-SI photoreceiver consists of an Si photodiode and a subsequent low-noise fixed gain amplifier. It is designed for ultra-fast conversion of small optical signals into equivalent output voltages. Operation is mostly self-explanatory. If in doubt, consult this document or contact support@femto.de.

For safe operation, please refer to the damage thresholds specified in the "Absolute Maximum Ratings", "Temperature Range" and "Power Supply" sections of this document.

The operating environment must be free of smoke, dust, grease, oil, condensing moisture, and other contaminants that could affect the operation or performance.

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY

Ultra High Speed Photoreceiver with Si-PIN Photodiode

Available Versions

HSA-X-S-1G4-SI-FST



1.035"-40 threaded flange with internally threaded coupler ring (outer diameter 30 mm) for free space applications. Compatible with many optical standard accessories and for use with various types of fiber connector adapters.

Optionally available:

Fiber adapters PRA-FC, PRA-FCA and PRA-FSMA. With the relative large 0.4 mm dia. photodiode installed in the HSA-X-S-1G4-SI input coupling is not critical. However, standard SM 9/125 fibers (PC or APC) with low numerical aperture (NA) are recommended for ensuring near 100% coupling efficiency.

HSA-X-S-1G4-SI-FC



Fix/permanent FC fiber connector for high coupling efficiency and excellent conversion gain accuracy.

Related Models

HSPR-X-I-1G4-SI-FST

Si-PIN, Ø 0.4 mm, 320 – 1000 nm, 1.4 GHz, inverting output, free space input, 1.035"-40 threaded flange

HSPR-X-I-1G4-SI-FC

Si-PIN, integrated ball lens, 320 – 1000 nm, 1.4 GHz, inverting output, FC fiber connector (fix/permanent)

HSA-X-S-2G-IN-FST

InGaAs-PIN, Ø 0.1 mm, 900 − 1700 nm, 2 GHz,

HSA-X-S-2G-IN-FC

IngaAs-PIN, ∅ 0.1 mm, 900 – 1700 nm, 2 GHz, free space input, 1.035"-40 threaded flange

InGaAs-PIN, integrated ball lens, 900 – 1700 nm, 2 GHz,

FC fiber connector (fix/permanent)

HSPR-X-I-2G-IN-FST

InGaAs-PIN,Ø 0.1 mm, 900 − 1700 nm, 2 GHz, inverting

output, free space input, 1.035"-40 threaded flange

HSPR-X-I-2G-IN-FC

InGaAs-PIN, integrated ball lens, 900 – 1700 nm, 2 GHz, inverting output, FC fiber connector (fix/permanent)

Available Accessories

PRA-FCA PRA-FSMA







Fiber-adapter with external 1.035"-40 thread (suitable for FST models only).

PS-15-25-L



Power Supply input: 100 – 240 VAC output: ±15 VDC

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY

E M T O

Page 2 of 6

Datasheet

HSA-X-S-1G4-SI

Ultra High Speed Photoreceiver with Si-PIN Photodiode

Specifications $V_S = +15 \text{ V}, T_A = 25 \text{ °C}, \text{ output load impedance } 50 \Omega,$

warm-up 20 minutes (min. 10 minutes recommended)

Gain Transimpedance gain 5.0×10^3 V/A (@ output load 50 Ω)

Conversion gain 2.55×10^3 V/W typ. (@ 760 nm, output load 50 Ω)

Frequency Response Lower cut-off frequency (–3 dB) 10 kHz

Upper cut-off frequency (–3 dB) 1.4 GHz (±15%)

Time Response Rise/fall time (10 % - 90 %) 250 ps (\pm 15%)

Input Noise equivalent power (NEP) 32 pW/√Hz (@ 760 nm, 100 MHz)

Optical saturation power 370 µW AC (for linear amplification, @ 760 nm) 10 mW CW (to prevent saturation, @ 760 nm)

Detector Si-PIN photodiode
Active area (FST version) Ø 400 μm

Active area (FC version) integrated ball lens

suitable for fibers up to 400 μm core diameter

Spectral range 320 – 1000 nm Max. sensitivity 0.51 A/W typ. (@ 760 nm)

Output Output voltage range $1.9 \text{ V peak-peak } (@ 50 \Omega \text{ output load})$

for linear operation and low harmonic distortion

 $\begin{array}{ll} \mbox{Output reflection S22} & -7.3 \mbox{ dB (@ f < 2.5 GHz)} \\ \mbox{Output impedance} & 50 \ \Omega \mbox{ (terminate with 50 } \Omega \mbox{ load)} \end{array}$

Output noise 3.6 mV RMS (24 mV peak-peak) typ. (@ 50 Ω load,

no signal on detector, measurement bandwidth 4 GHz)

Optical Input Connector Material FST flange 1.4305 stainless steel, nickel-plated

Material FST coupler ring 1.4305 stainless steel, glass bead blasted

Material FC receptacle nickel silver

Power Supply Supply voltage +15 V

Supply current 130 mA (depends on operating conditions,

recommended power supply capability min. 200 mA)

Case Weight 133 g (0.29 lbs) HSA-X-S-1G4-SI-FST incl. coupler ring

110 g (0.24 lbs) HSA-X-S-1G4-SI-FC

Material AlMg4.5Mn, nickel-plated

Temperature Range Storage temperature -30 °C ... +85 °C

Operating temperature 0 °C ... +60 °C

Absolute Maximum Ratings Optical input power (CW) 12 mW (averaged)

Power supply voltage 20 V

OVERY F E M T O

Datasheet

HSA-X-S-1G4-SI

Ultra High Speed Photoreceiver with Si-PIN Photodiode

Connectors Input HSA-X-S-1G4-SI-FST 1.035"-40 threaded flange for

free space applications and for use with various types of optical

standard accessories


HSA-X-S-1G4-SI-FC FC fiber optic connector

(fix/permanent, FC/PC and FC/APC compatible)

Output SMA jack (female)

Power supply LEMO® series 1S, 3-pin fixed socket

(mating plug type: FFA.1S.303.CLAC52)

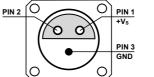

Pin 1: +15 V

Pin 2: NC Pin 3: GND

HSA-X-S-1G4-SI, internally threaded coupler ring (FST version only), LEMO® 3-pin connector, datasheet, transport package

HSA-X-S-1G4-SI-FST 1.035"-40 threaded flange for free space applications and

for use with various types of optical standard accessories.


HSA-X-S-1G4-SI-FC FC fiber optic connector

(fix/permanent, FC/PC and FC/APC compatible).

Spectral Response

Scope of Delivery

Ordering Information

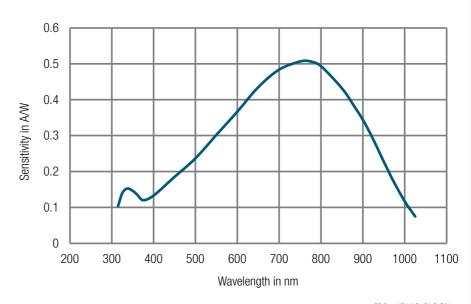


DB-Sens-HSA-X-S-1G4-SI_R01

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY

Ultra High Speed Photoreceiver with Si-PIN Photodiode

Dimensions

HSA-X-S-1G4-SI-FST (1.035"-40 threaded free space input)

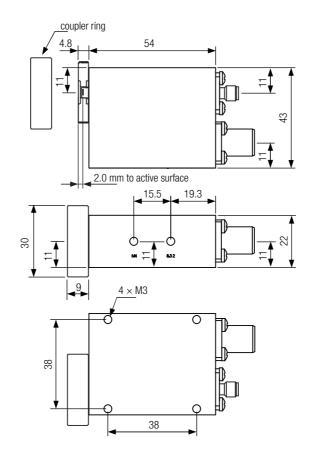

DZ-HS-FST_R1


all dimensions in mm unless otherwise noted


SOPHISTICATED TOOLS FOR SIGNAL RECOVERY


Ultra High Speed Photoreceiver with Si-PIN Photodiode

Dimensions (continued)

HSA-X-S-1G4-SI-FC (FC fiber optic connector)

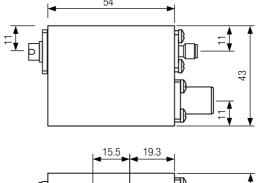

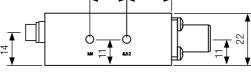

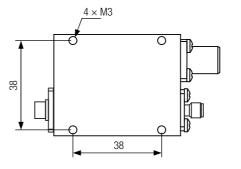

DZ-HS_FC_R1

all dimensions in mm unless otherwise noted

FEMTO Messtechnik GmbH Klosterstr. 64 10179 Berlin · Germany Phone: +49 30 280 4711-0 Fax: +49 30 280 4711-11 Email: info@femto.de www.femto.de Specifications are subject to change without notice. Information provided herein is believed to be accurate and reliable. However, no responsibility is assumed by FEMTO Messtechnik GmbH for its use, nor for any infringement of patents or other rights of third parties which may result from its use. No license is granted by implication or otherwise under any patent or patent rights of FEMTO Messtechnik GmbH. Product names mentioned may also be trademarks used here for identification purposes only.

© by FEMTO Messtechnik GmbH · Printed in Germany

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY